


GoldiBlot™

See His-tagged proteins in ~1 hour
with Ni-NTA-Nanogold® Technology

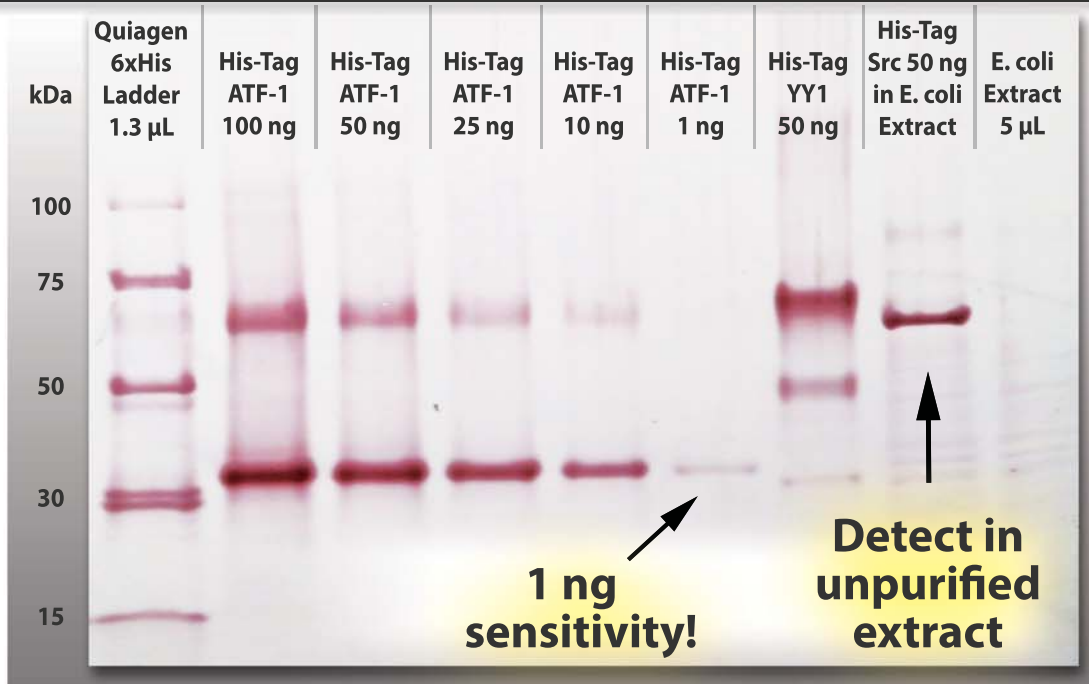

Ni-NTA binds directly to His-tags *No antibodies needed!*

- **Fast** Results in about an hour!
- **Directly visualized by eye**
No film, no equipment required!
- **Colorimetric detection**
Low background - Dark, clean signal
- Highly **Sensitive** ...to 1 ng!
- **Saves money**
No more expensive antibodies!

No need to purify...
Just load crude extract!

Transferred membrane	
1. TBST wash	3 min.
2. Block membrane	15 min.
3. GoldiBlot™ Incubation	30 min.
4. Imidazole-TBST wash	2 x 3 min.
5. dH ₂ O wash	3 x 3 min.
6. GoldiBlot™ AutoMet detection	10 min.

TOTAL TIME: 1 HOUR 13 min.

Available in
5-blot trial size
or 15-blot kits



Tel: 631-205-9490 Fax: 631-205-9493
US toll free: 877-447-6266

Questions? Ask our scientists: tech@nanoprobes.com